

Tuning the Collimating lens.

When you receive your collimator it will be set up for your microscope. For fine tuning please follow the instructions below.

Close shutter on Lumen 200.

Loosen the liquid light guide locking screw and turn part A to focus beam, lock with the locking screw.

Open shutter and check beam alignment.

Repeat as necessary.

DO NOT COMPLETE THIS WITH LIGHT EMITTING FROM THE LIGHT GUIDE, WARNING HIGH LEVELS OF LIGHT ARE EMITTED FROM THE LUMEN 200 LIGHT GUIDE.

Figure a. Microscope adaptor and collimator.

